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Safety Instructions Lab Burner „Fireboy Plus“ (Integra)

1. Before handling of the Fireboy Plus you have to make sure that the black screw on top of the burner connected cartridge adapter (red arrow in photo below, burner unit and cartridge not visible) is in the “ON” position completely, that means tightened CCW. *This screw is only allowed to be turned when changing cartridges.*

2. When handling the Fireboy Plus you have to keep a *safety distance* of app. half of an arm's length. You must wear safety glasses. Generally there is the risk of burn because of the open flame and hot surfaces.

3. The burner can be switched on by pressing the „menu“ button for ap. 2 sec. Display gets illuminated and the word “Button” appears. Pressing the menu button once again shortly starts the burner. Height and temperature of the flame can be adjusted by the two buttons left and in the middle. Press again menu button shortly to switch flame off. *After finishing your work you must switch off the burner completely by pressing the menu button for app. 2 sec (Display goes out).*

For safety precautions the Fireboy Plus *must* be used in the “Button” mode only. If a different word appears in the display turn menu button until “Button” appears.

4. *Transport of burner in the lab:* For safety reasons the burner is only allowed to be moved in a cold and completely switched off state (display completely off).

5. *Exchanging the CV360 cartridge:* Fireboy Plus must be cold and switched off completely (Display off). Black screw from cartridge adapter can be turned CW in the “OFF” position (cartridge exchange position) than. Remove empty cartridge from adapter by pulling firmly. Put in new cartridge into adapter again by overcoming a drag, hear a “click” again. Turn the black screw of the adapter in the “ON” position (working mode) CCW again. The cartridge with the open valve is now connected to the burner again.

6. *Cleaning burner:* Fireboy Plus should be cleaned on a frequent basis according to the manual.

In case of any problems please contact your group leader, Jan Brix or Rainer Liefländer.